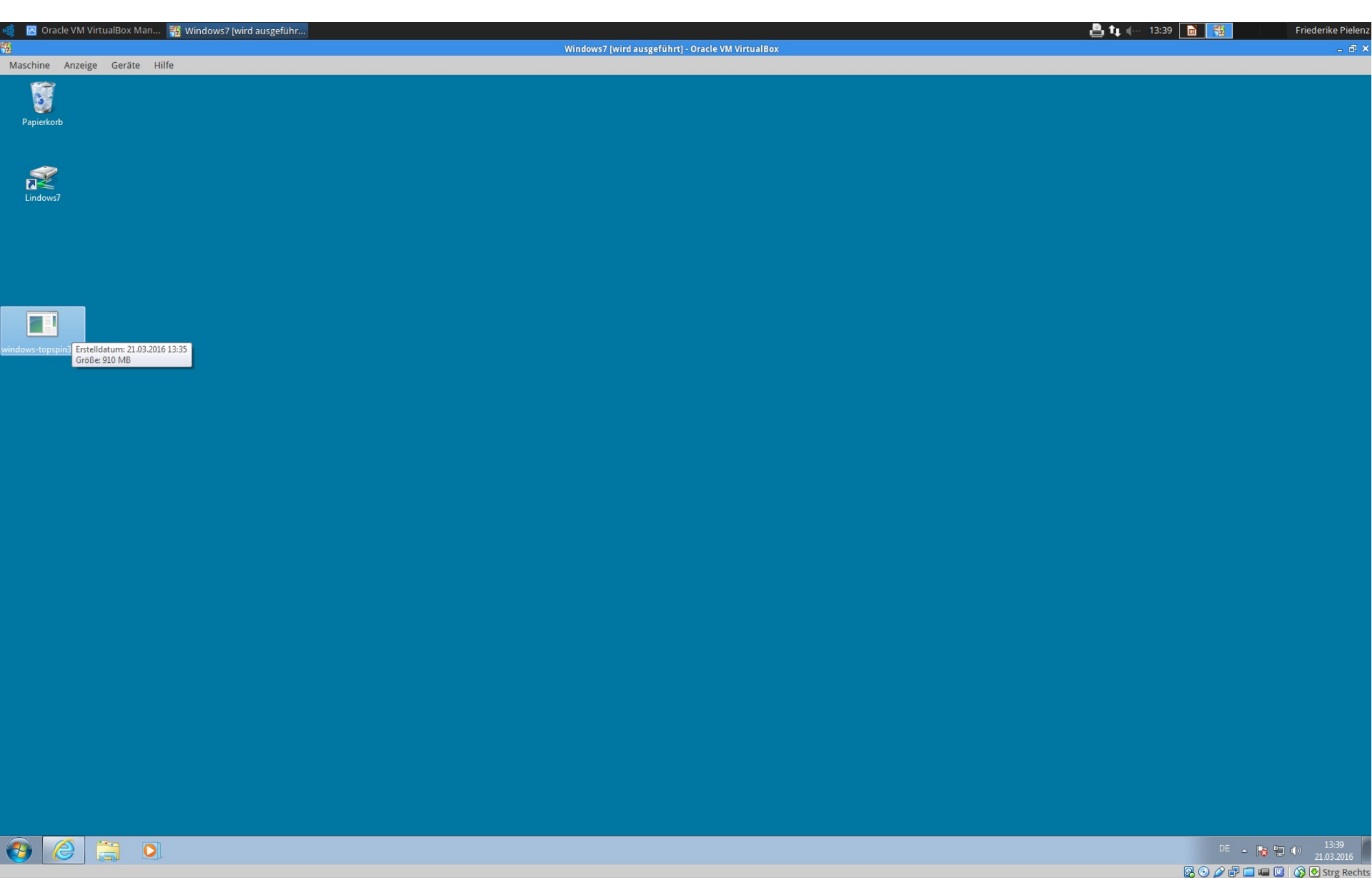
Installationsanleitung TopSpin (Windows 7)

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install@EDV

BRUKER

Bruker NMR Software Installer
including TopSpin 3.5pl2 and CMC-a

Overall progress

- ✓ Setup parameters complete
- ✓ Check consistency complete
- Select target directories running
- Prepare harddisk -
- Install files -
- Post processing -

Status: Selecting installation directory for TopSpin

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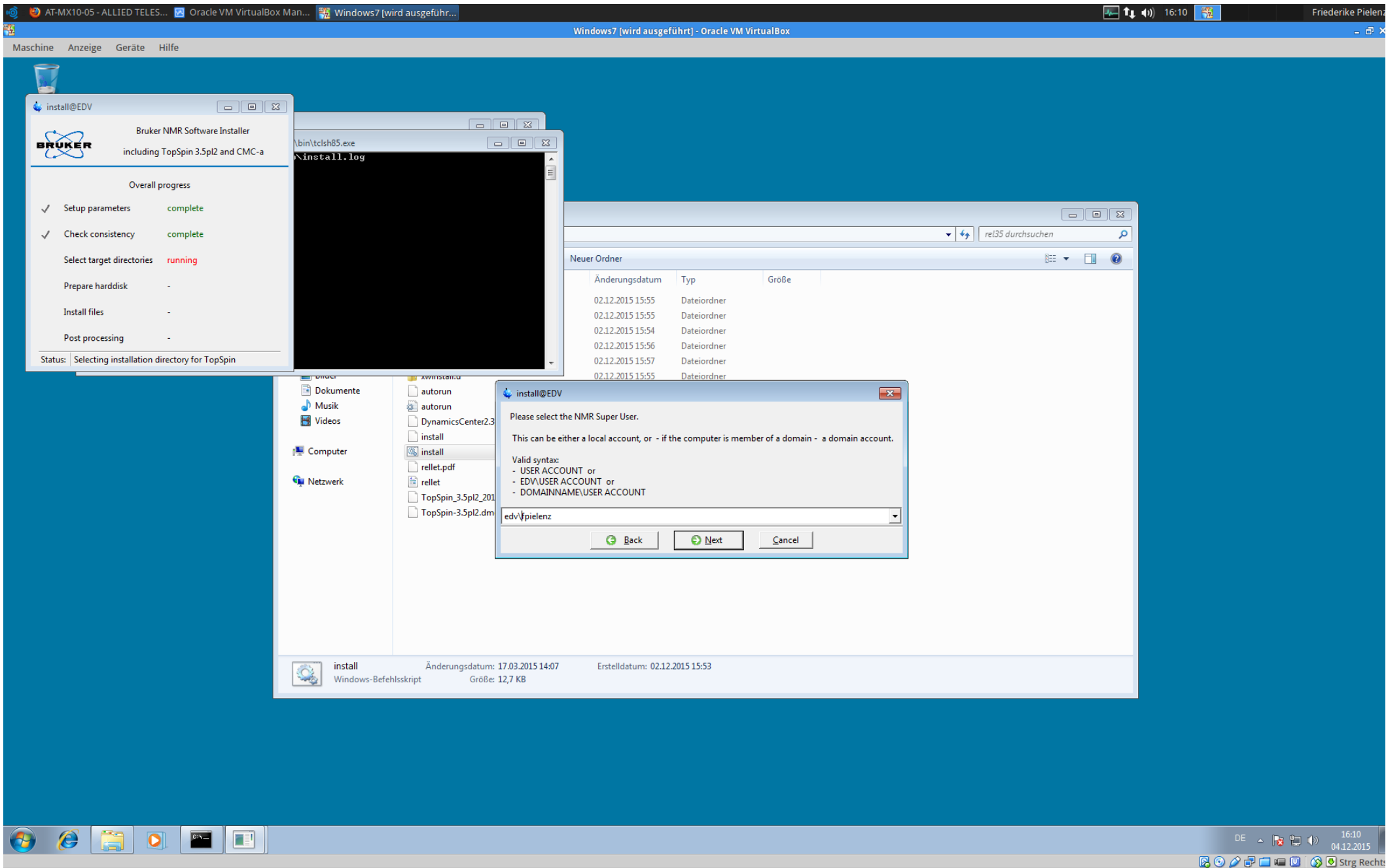
install@EDV

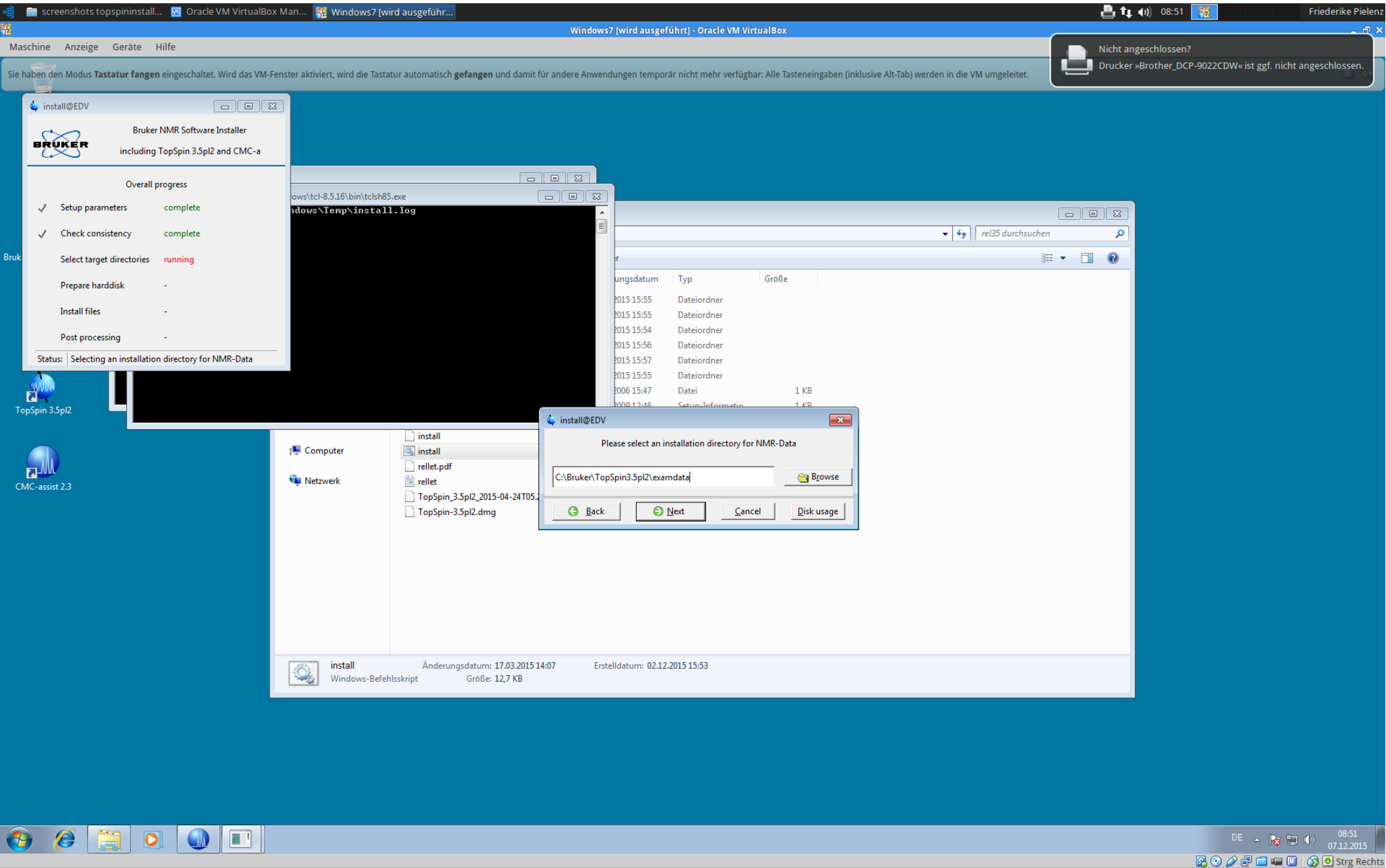
Please choose an NMR Administration Password
(requested for commands like 'cf', 'expinstall' etc.)

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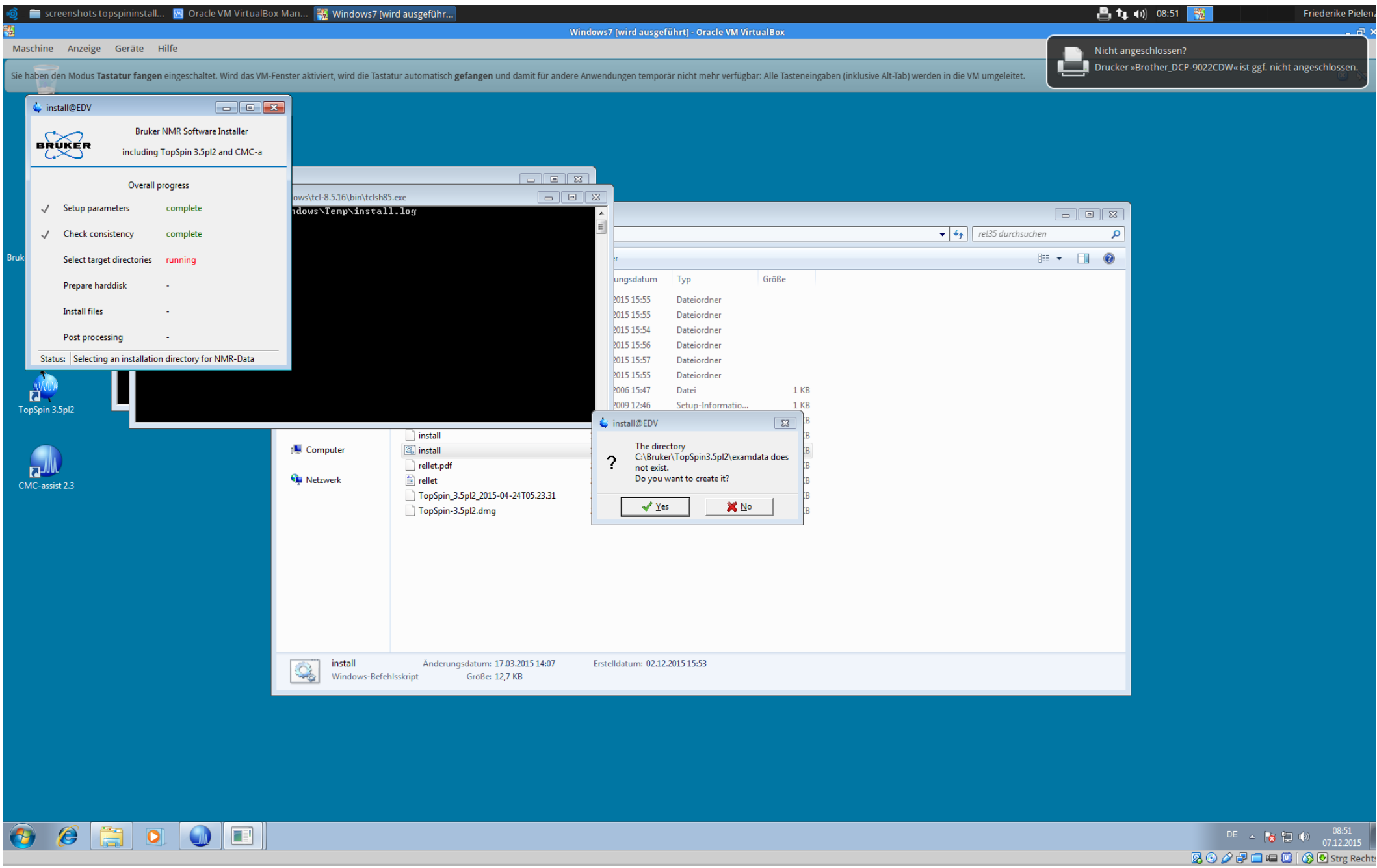
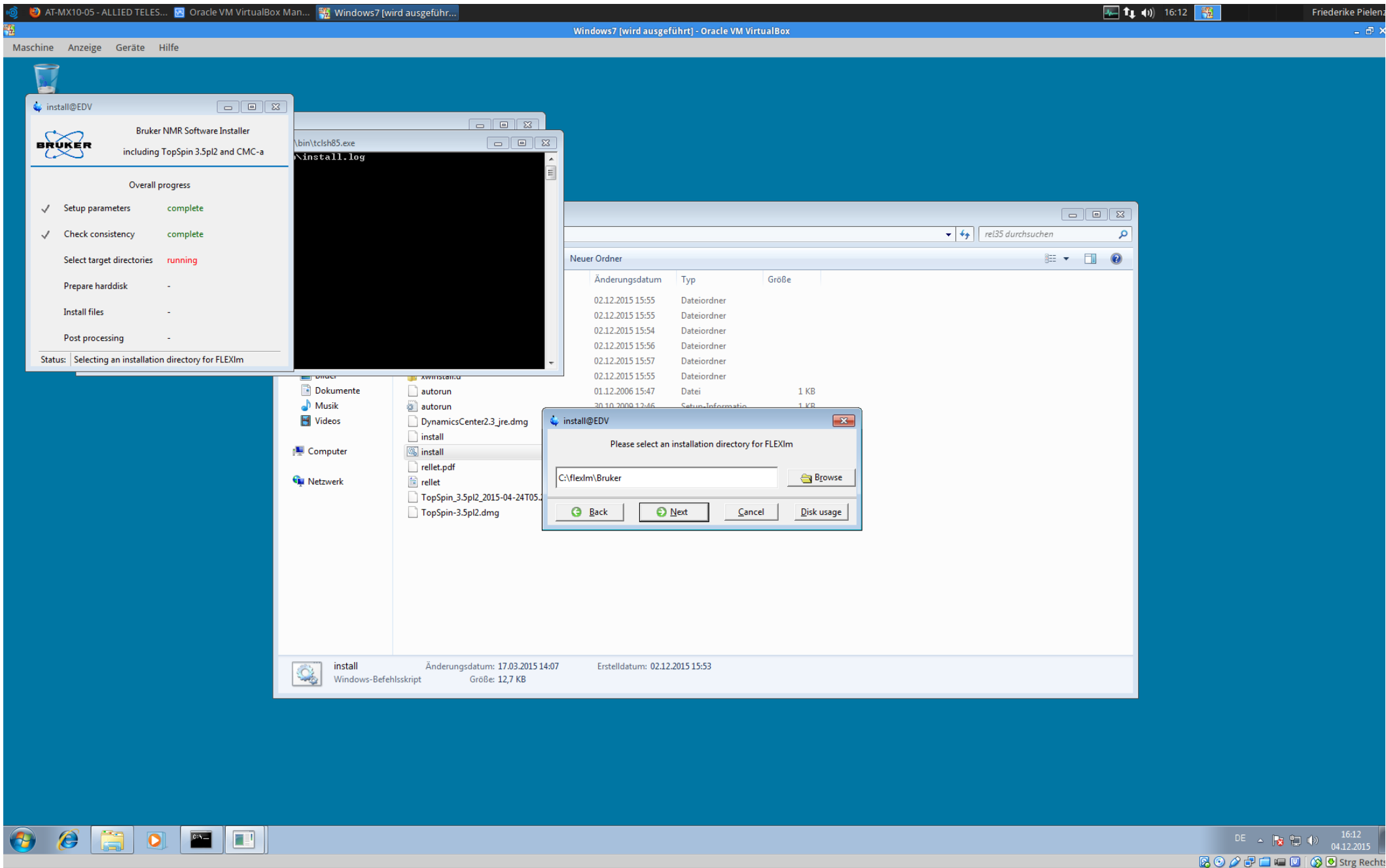
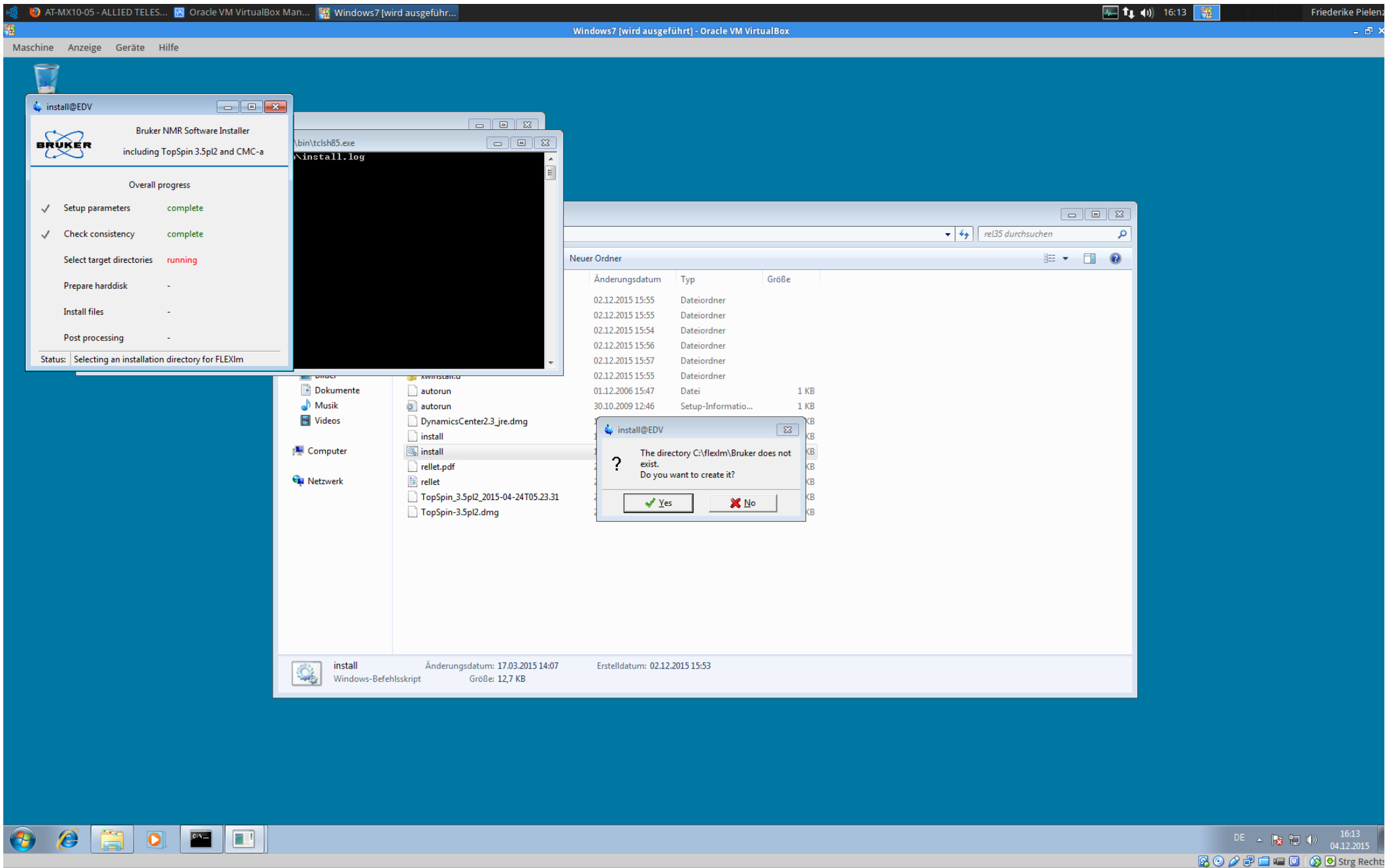
install@EDV

Windows firewall is installed on your system.

If it has not been disabled, starting FLEXIm and TopSpin will bring up several information dialogs about 'blocked components'. To prevent this, the firewall must be configured to allow FLEXIm and TopSpin operations.

The installation can do this now, or an Administrator can do it any time later, running

```
C:\Windows\system32\cmd.exe /c C:\Bruker\TopSpin3.5pl2\prog\bin\Utilities\Miscellaneous\setfirewall.cmd install program
```

```
C:\Windows\system32\cmd.exe /c C:\Users\fpielanz\Desktop\rel35\windows\bin\setfirewall.cmd install flexim
```

Shall the installation configure the firewall for you?

Yes, configure system firewall automatically

[Automatically creates all firewall rules necessary for the products selected for installation](#)

Show necessary firewall rules


This is useful if you intend to change custom firewall rules manually.
The firewall rules will be shown after the products have been installed onto the hard disk.

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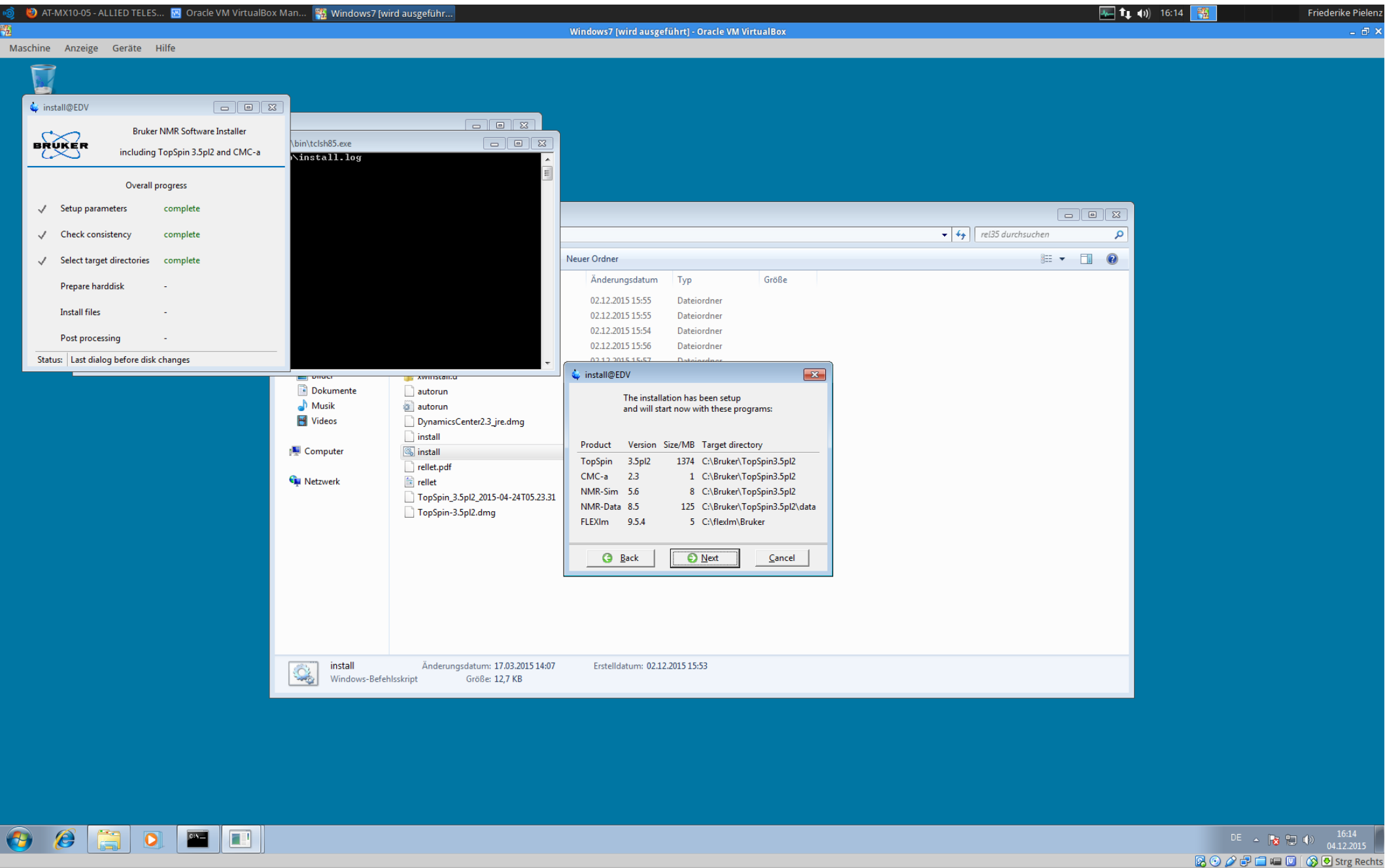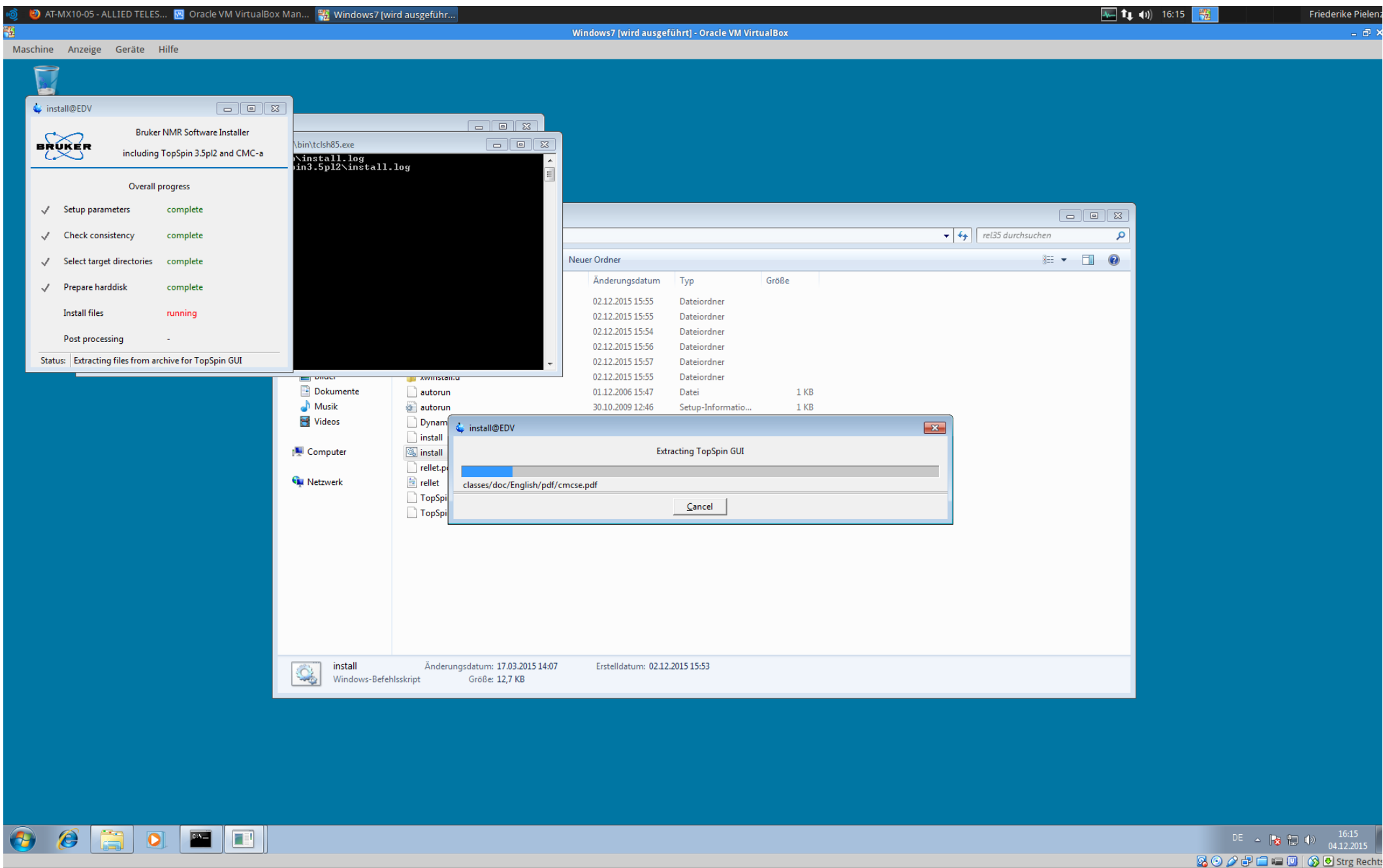
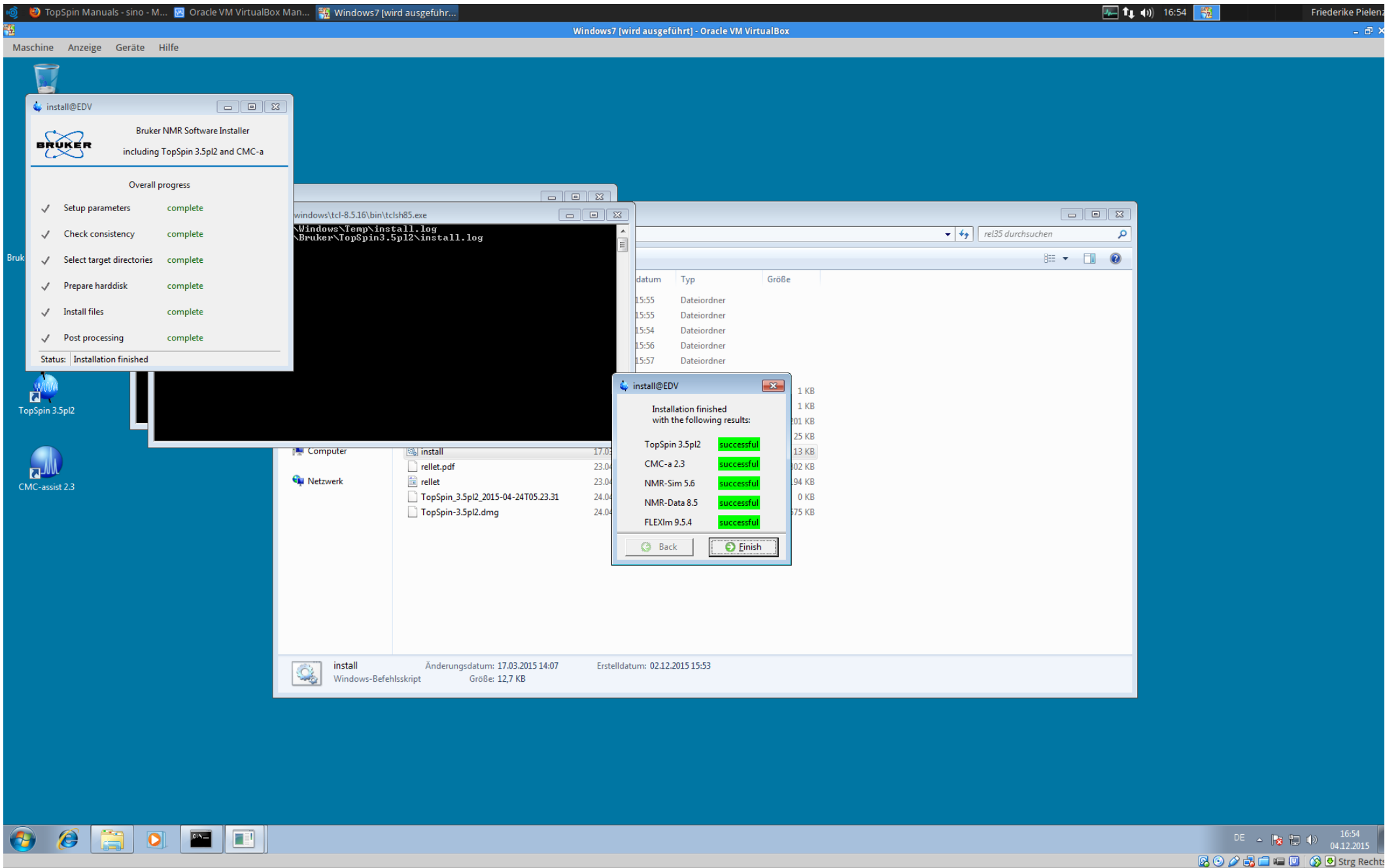
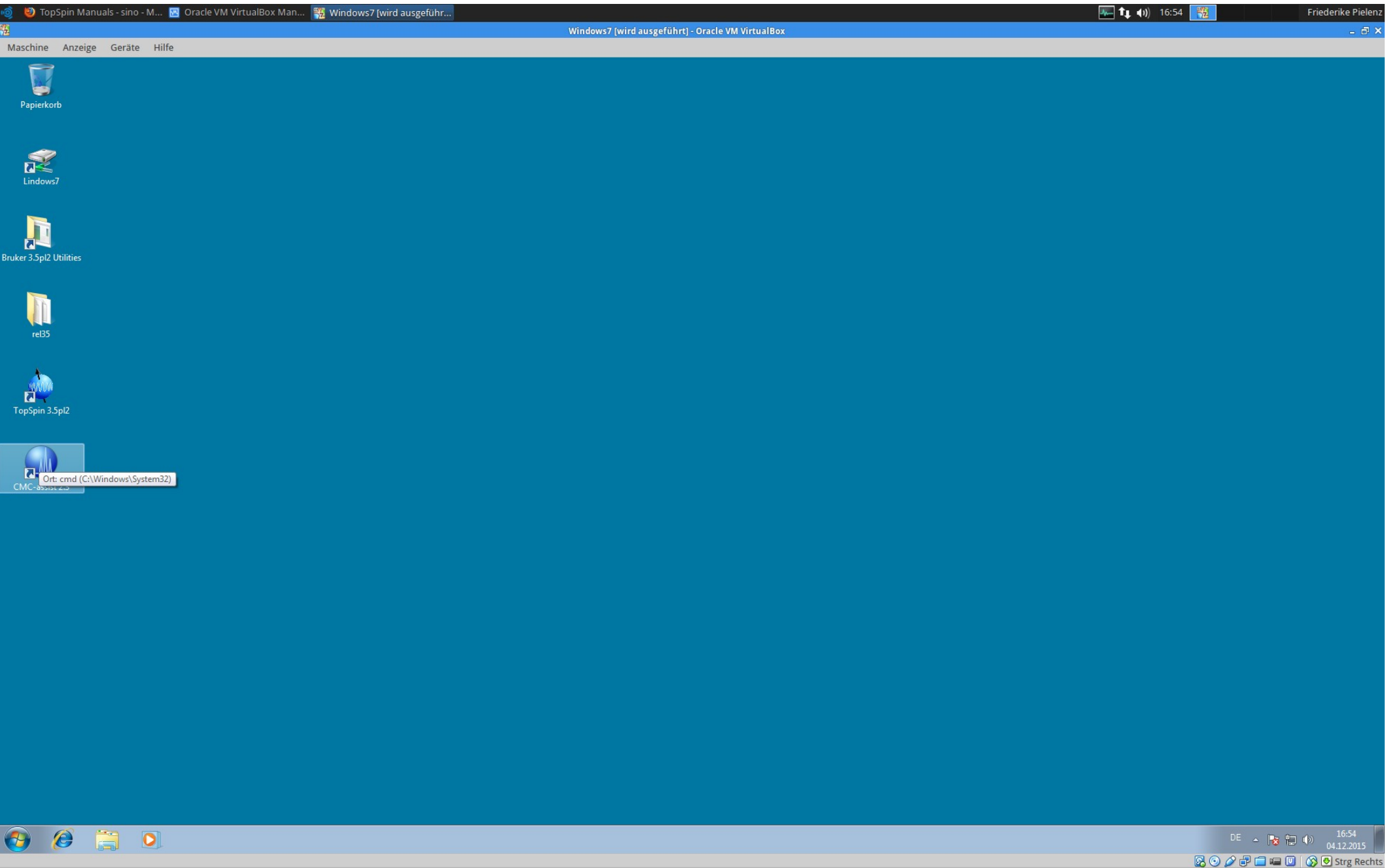
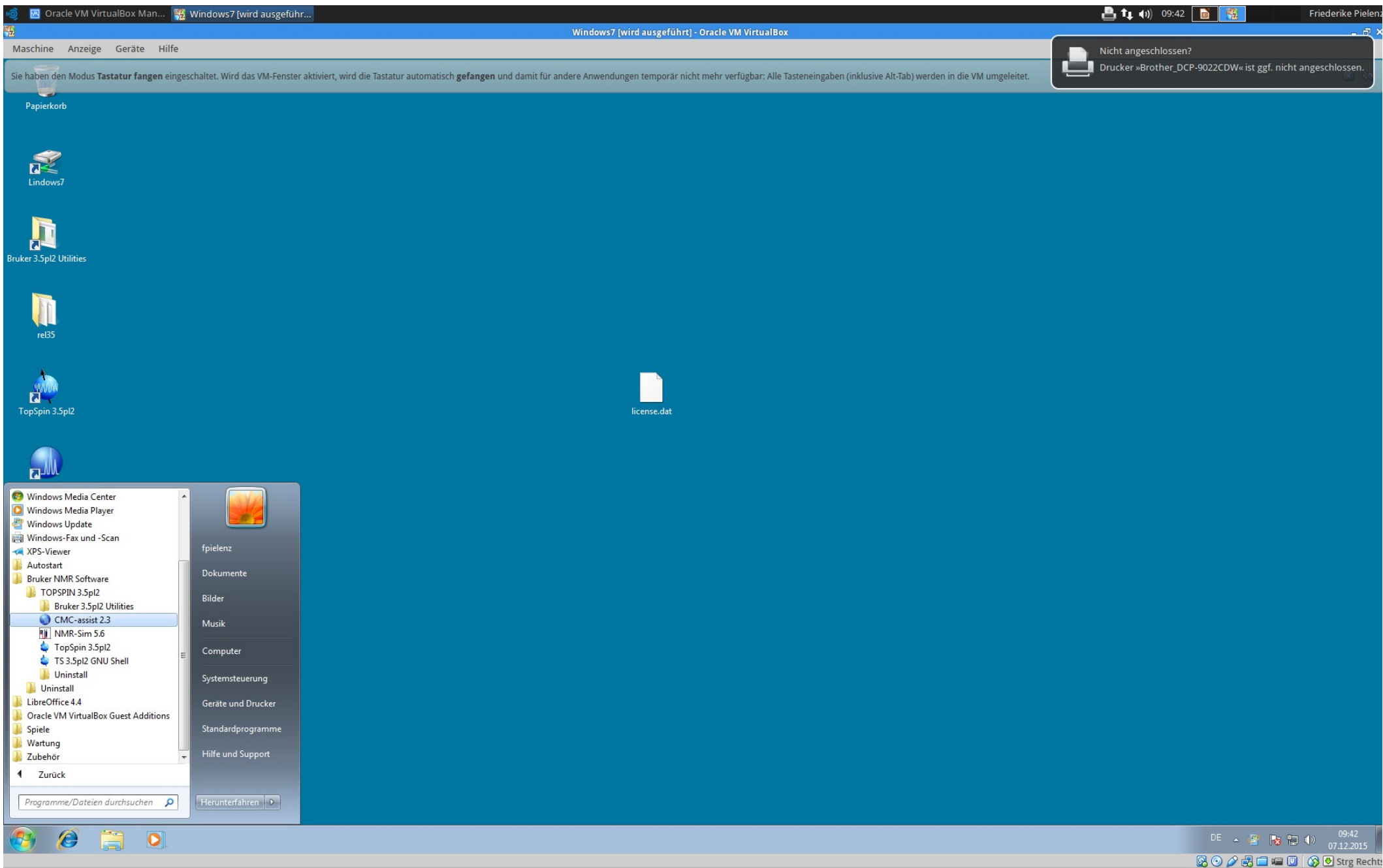
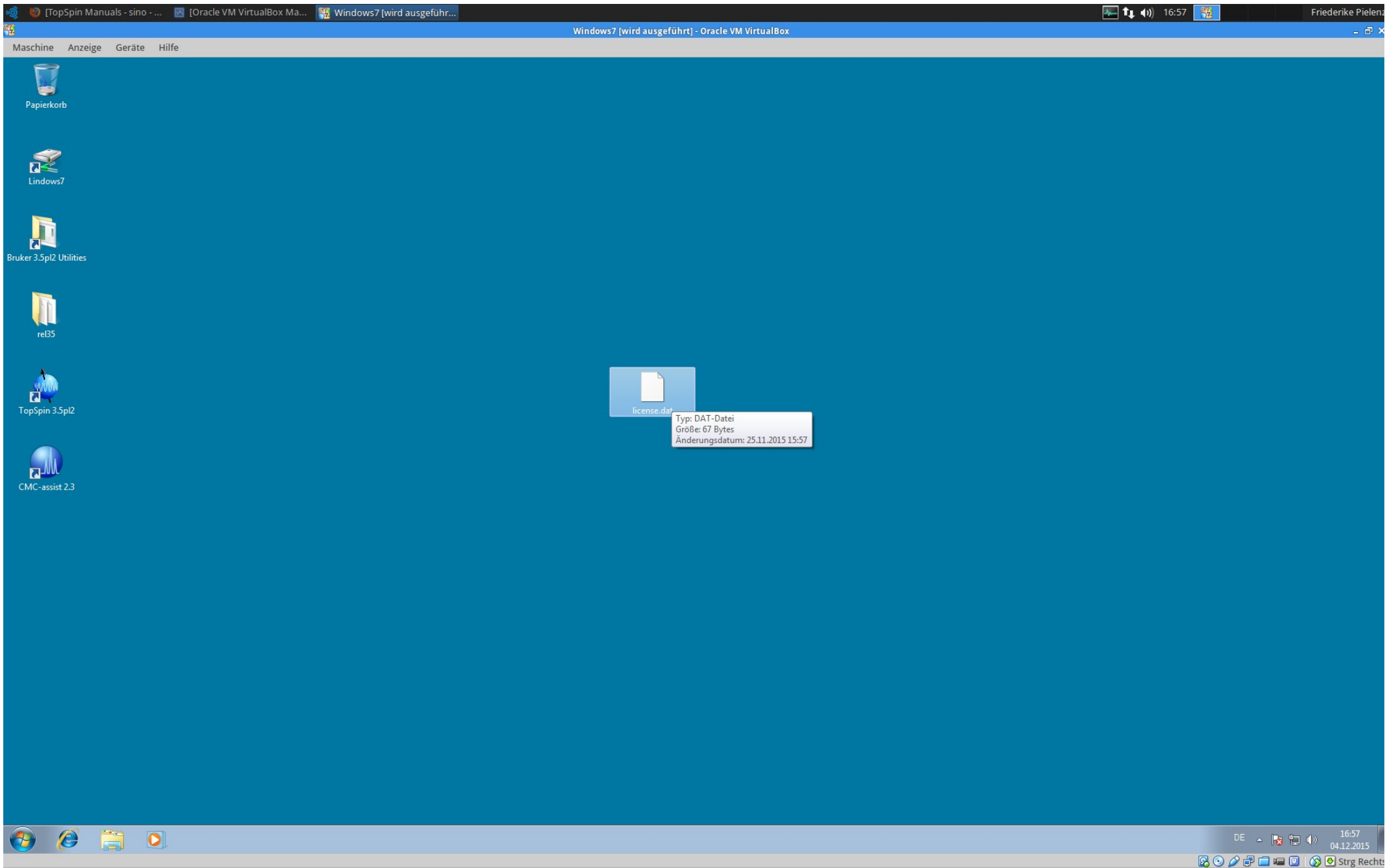
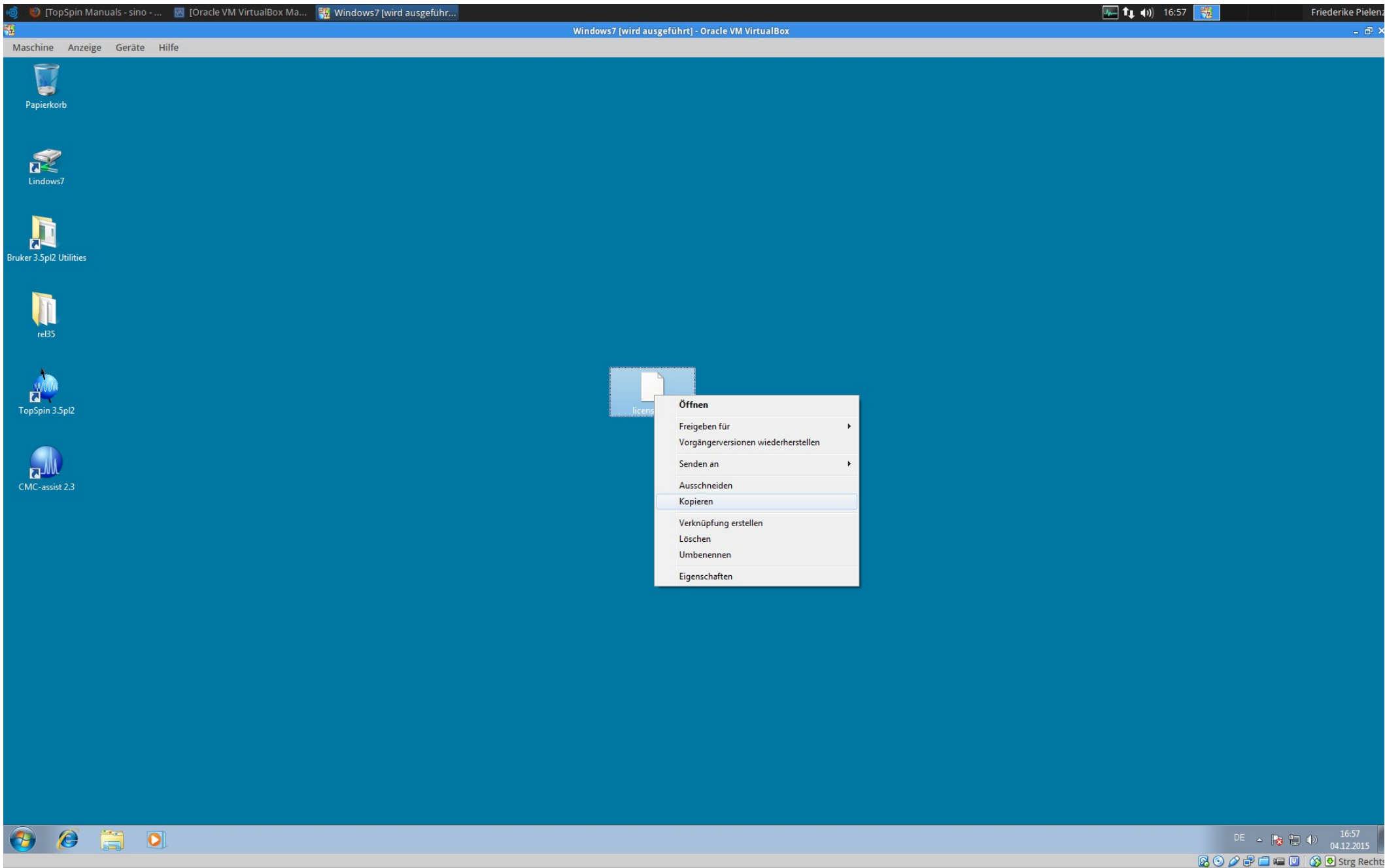
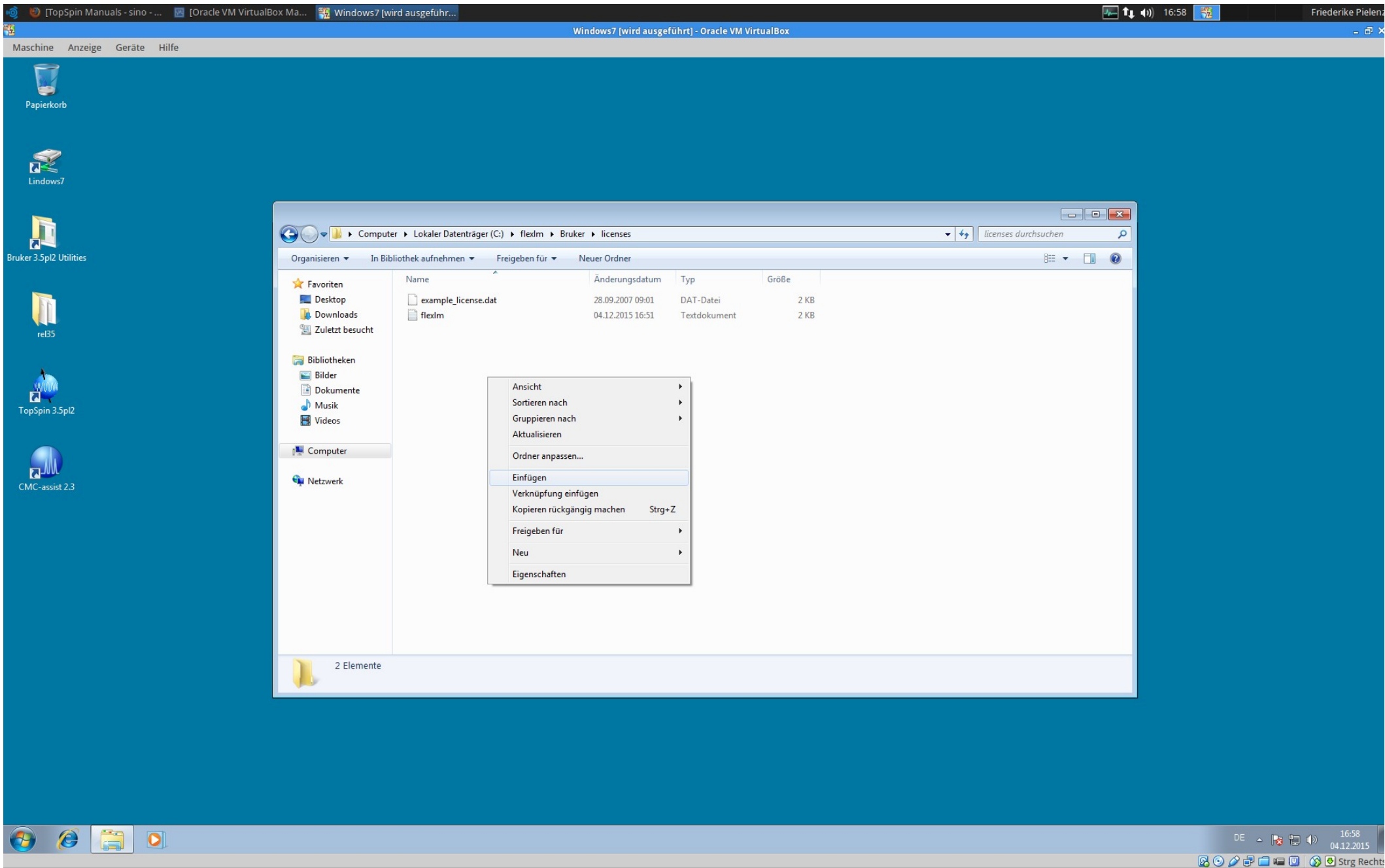
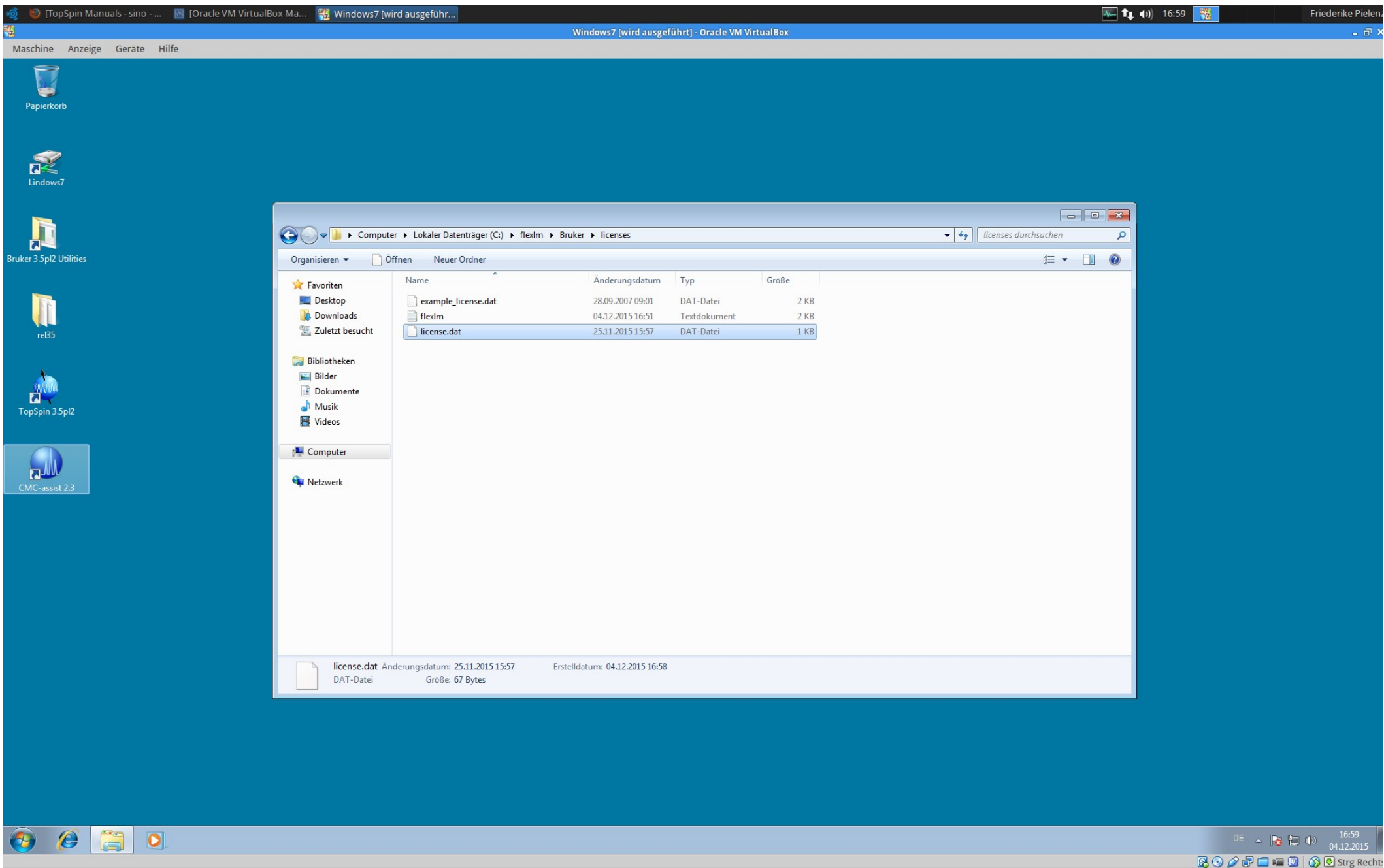
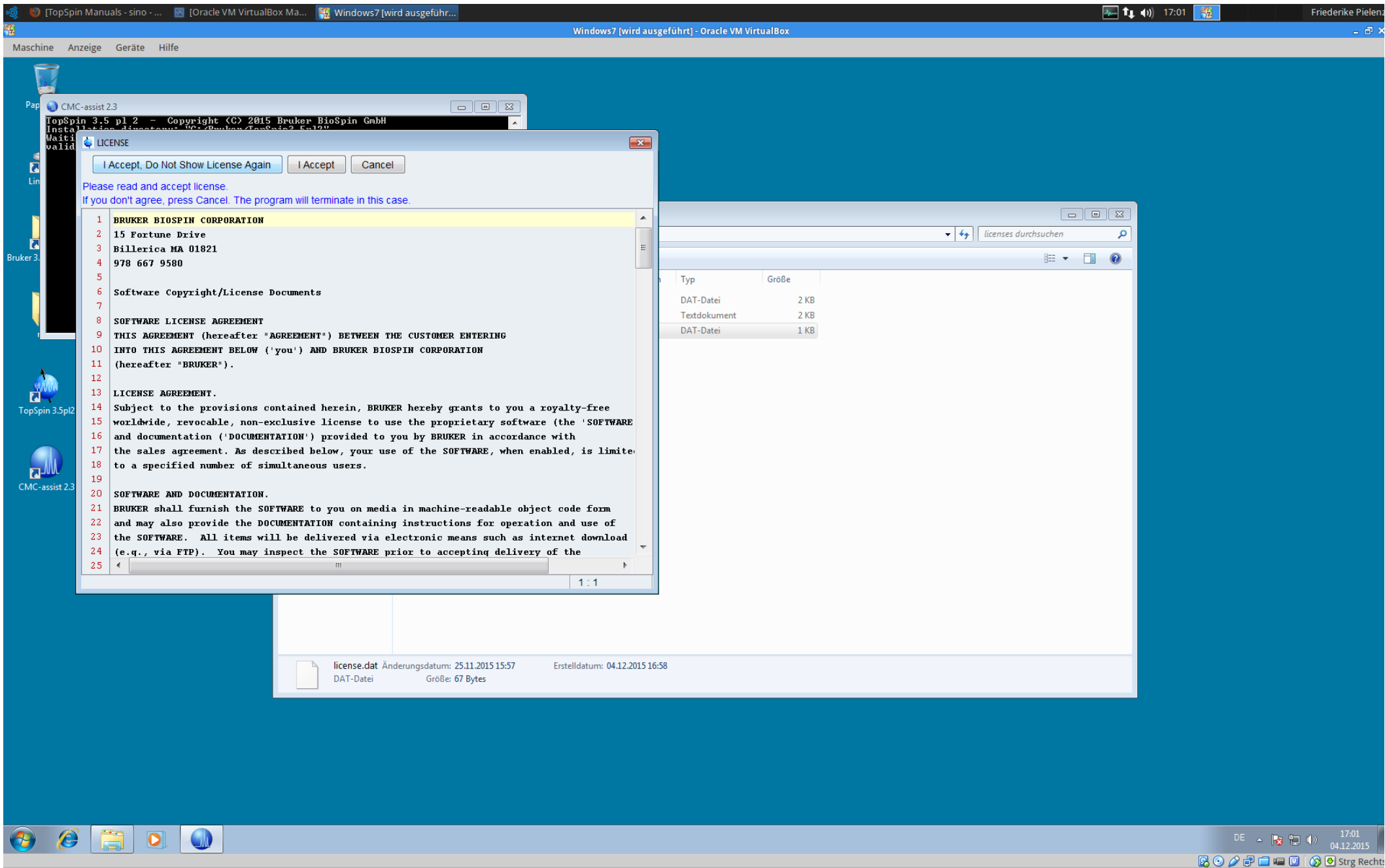
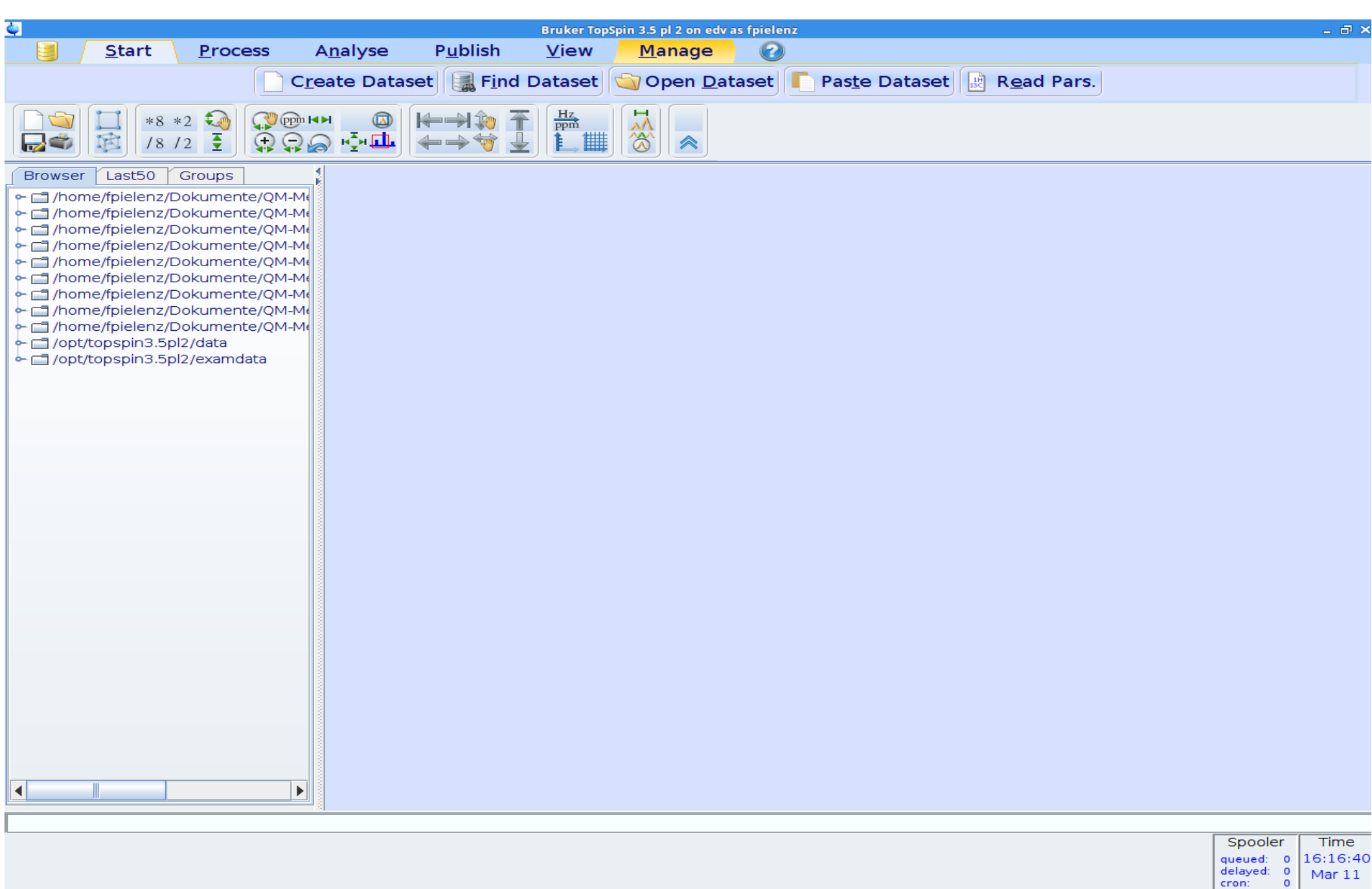
Bruker TopSpin 3.5 pl 2 on edv as fpielenz

Start Process Analyse Publish View **Manage** ?

Preferences Spectrometer Security Commands Remote


Setup user preferences
Allows you to change colors, fonts and many other properties influencing the behaviour of TopSpin.

Browser Last50 Groups

- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /home/fpielenz/Dokumente/QM-Me
- /opt/topspin3.5pl2/data
- /opt/topspin3.5pl2/examdata


Spooler	Time
queued: 0	16:17:44
delayed: 0	Mar 11
cron: 0	

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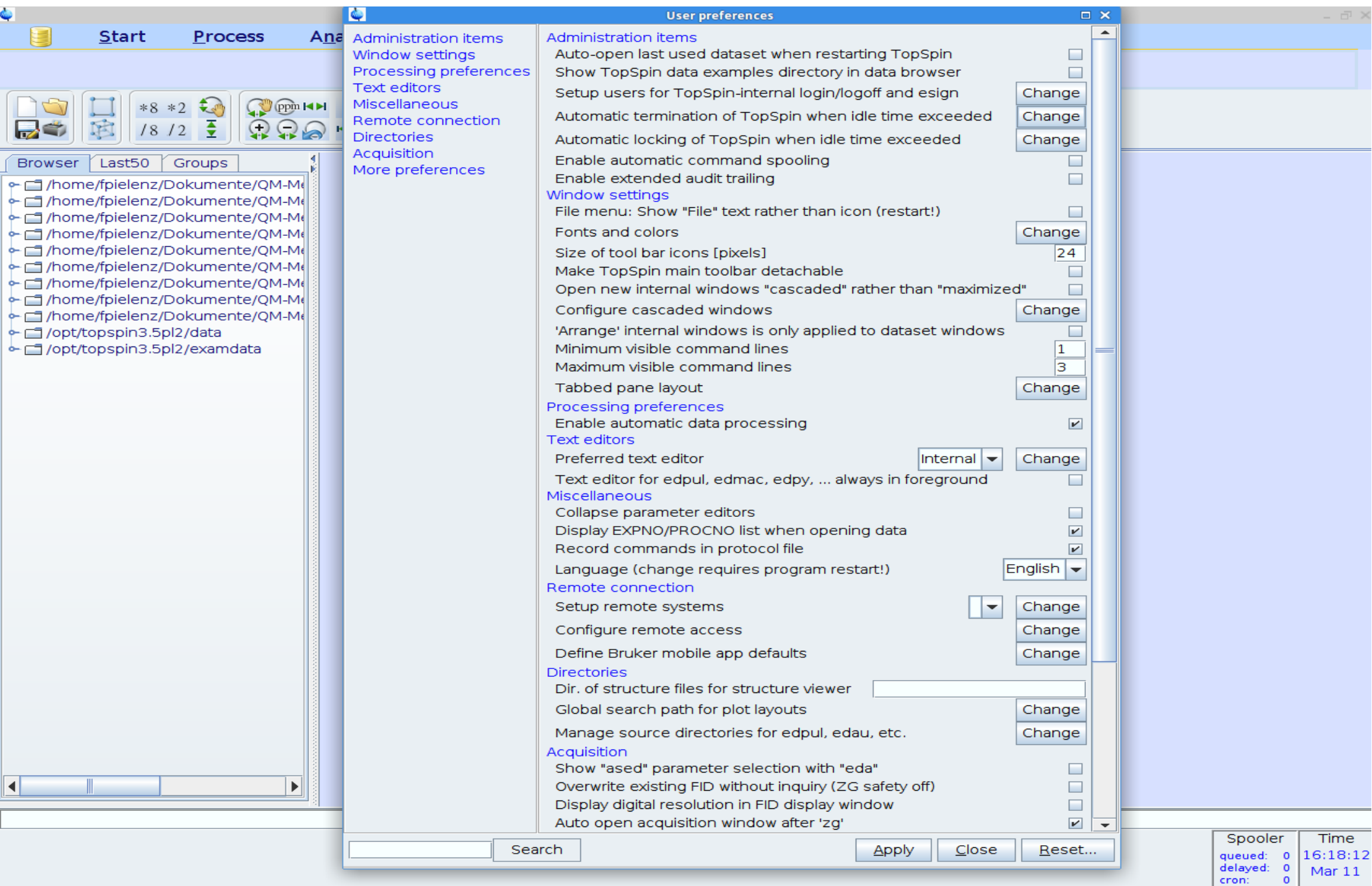Spooler	Time
queued: 0	16:18:12
delayed: 0	Mar 11
cron: 0	

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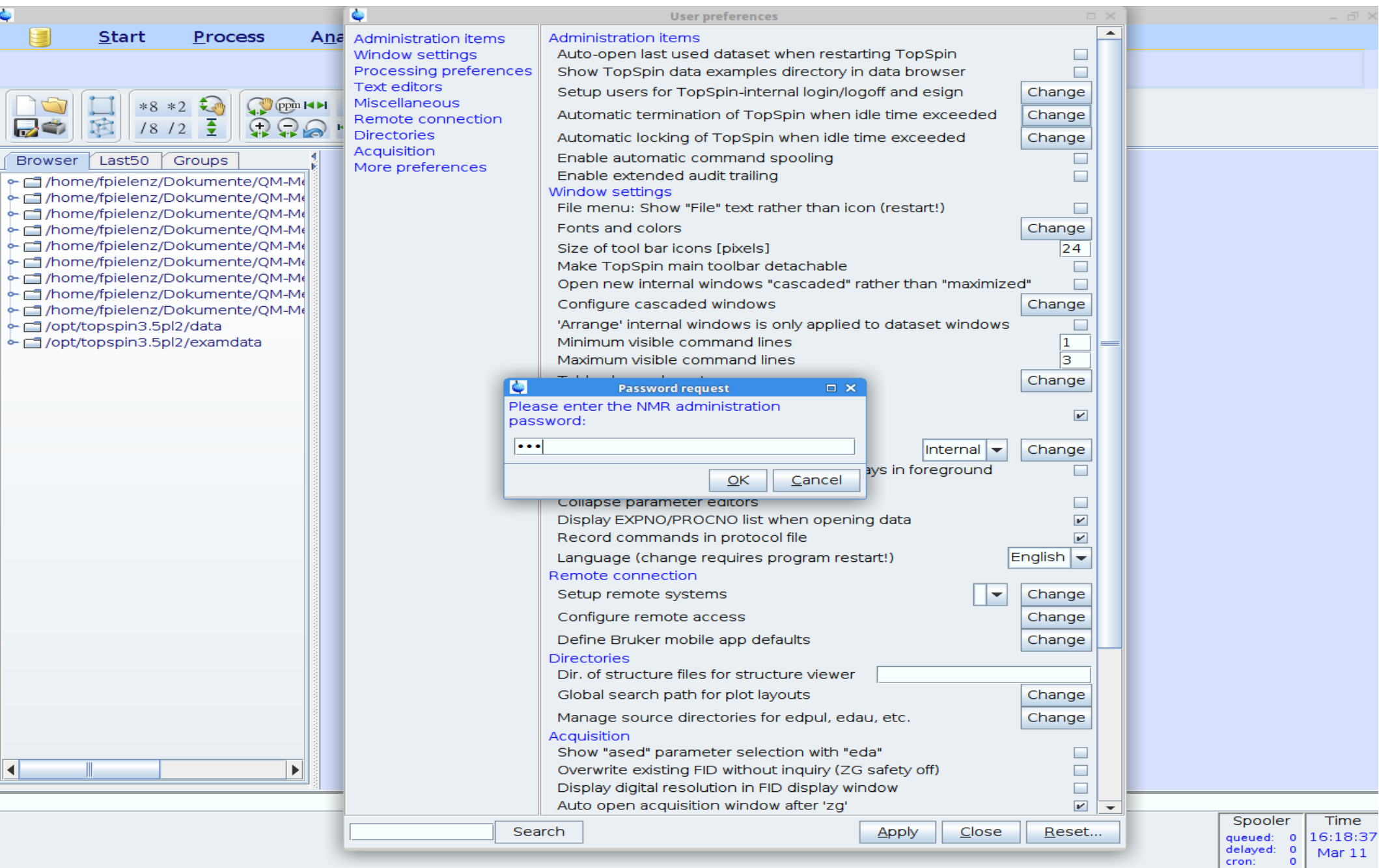The screenshot shows the 'User preferences' dialog box in the TopSpin software. A sub-dialog box is open, prompting the user to enter the maximum allowed idle time in minutes. The value '30' is entered in the input field. The background shows the 'Administration items' section of the preferences, including 'Automatic termination of TopSpin when idle time exceeded'.

Please enter the maximum allowed idle time in minutes. TopSpin will terminate when in idle state longer than this time. Enter zero to disable auto-termination. You must restart TopSpin to make changes effective.

30

OK Cancel

Administration items

- Auto-open last used dataset when restarting TopSpin
- Show TopSpin data examples directory in data browser
- Setup users for TopSpin-internal login/logoff and esign
- Automatic termination of TopSpin when idle time exceeded
- Automatic locking of TopSpin when idle time exceeded
- Enable automatic command spooling
- Enable extended audit trailing

Window settings

- File menu: Show "File" text rather than icon (restart!)
- Fonts and colors
- Size of tool bar icons [pixels] 24
- Make TopSpin main toolbar detachable
- Open new internal windows "cascaded" rather than "maximized"
- Configure cascaded windows
- 'Arrange' internal windows is only applied to dataset windows
- Minimum visible command lines 1
- Maximum visible command lines 3

Internal

background - Display EXPNO/PROCNO list when opening data
- Record commands in protocol file
- Language (change requires program restart!) English

Remote connection

- Setup remote systems
- Configure remote access
- Define Bruker mobile app defaults

Directories

- Dir. of structure files for structure viewer
- Global search path for plot layouts
- Manage source directories for edpul, edau, etc.

Acquisition

- Show "ased" parameter selection with "eda"
- Overwrite existing FID without inquiry (ZG safety off)
- Display digital resolution in FID display window
- Auto open acquisition window after 'zg'

Search Apply Close Reset...

Spooler	Time
queued: 0	16:19:05
delayed: 0	Mar 11
cron: 0	

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